

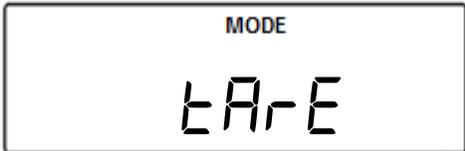
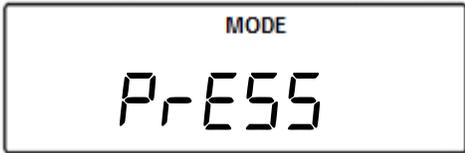
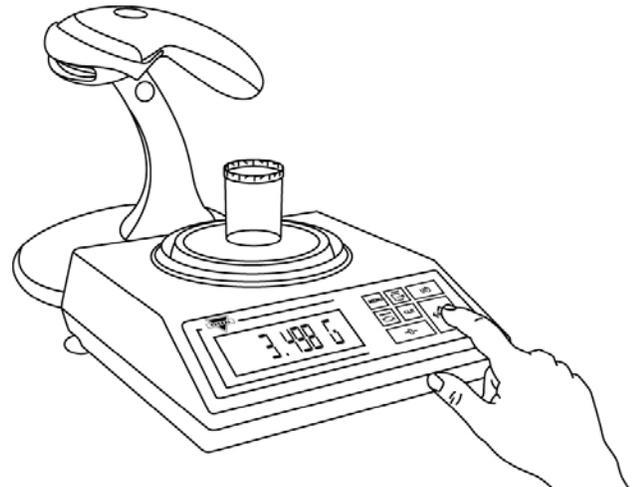


Counting Prescription Drugs for the first time (DRX-4C)



1. "P-r-E-S-S T-A-R-E" will appear on the scale's display, indicating to tare a counting container.

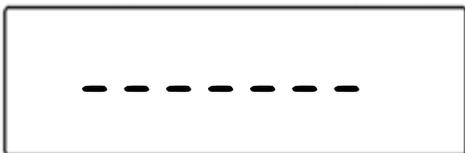


2. Place an empty vial or counting container on the pan. The container's weight will be displayed. Wait for the weight to stabilize and press the T key to tare.

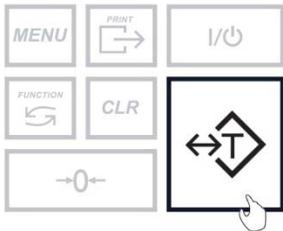



3. Once the scale has finished taring "S-c-A-n b-o-t-t-l-E" will appear on the display indicating you should scan the NDC code associated with the drug.

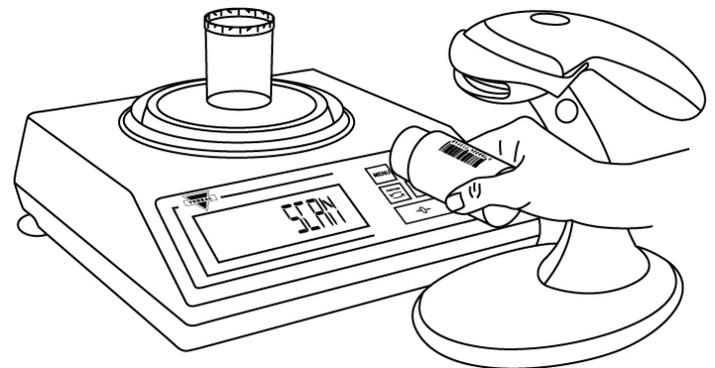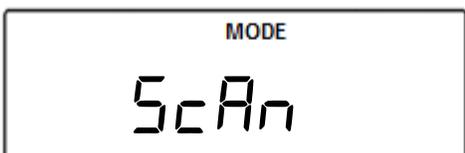

- Scan the NDC code by slowly placing it under the scanner light.
- The scale will display SAPL On (SAPL On). Manually count the required piece sample and place it in the container on the pan.

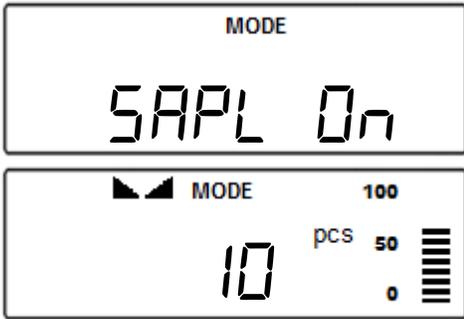


- The scale will display the weight of the sample. Press the T key to confirm the sample weight.

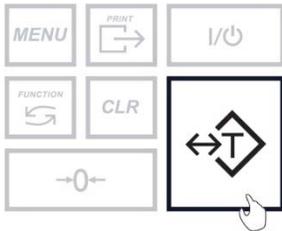
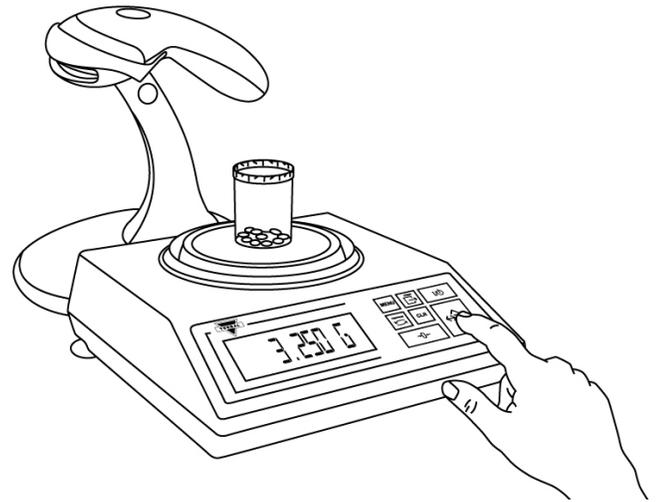


- The scale will display FILL indicating to begin filling the remainder of the count. Begin to pour the pills into the container.

- Once an accurate piece count has been taken, the container and its content may be removed from the scale. After removing the container, done will appear on the screen.

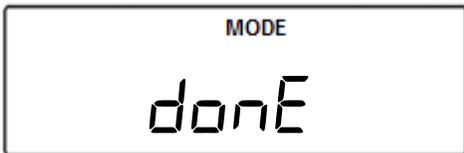


- To finish the count and begin a new one, press the Clear key.

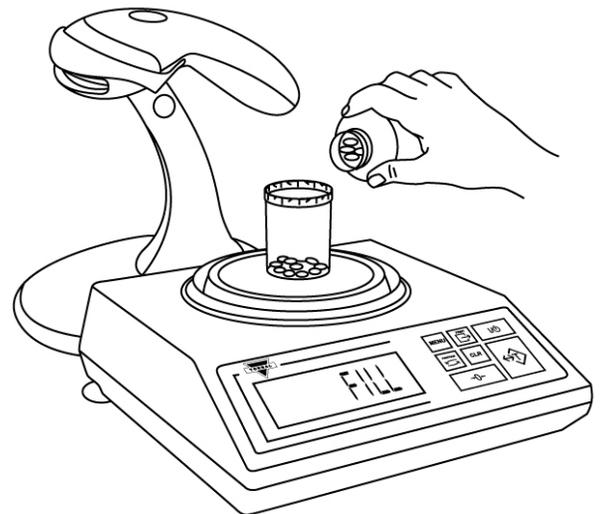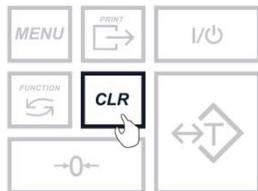


Counting Prescription Drugs Stored in the Database (DRX-4C Only)




- To enter Pill Counting Mode, press the function key. Select “count” by pressing the T key.

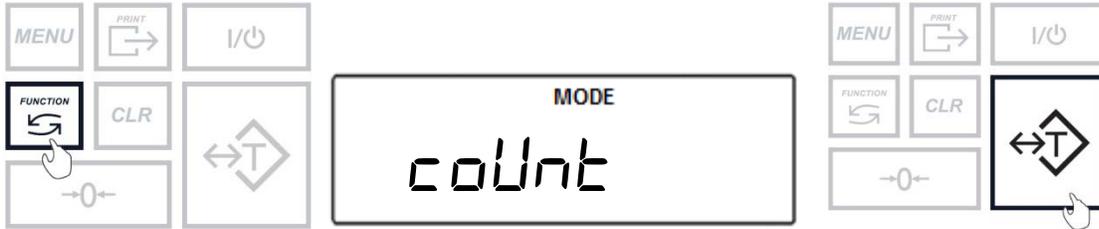



- “PRESS TARE” will appear on the scale’s display, indicating to tare a counting container.

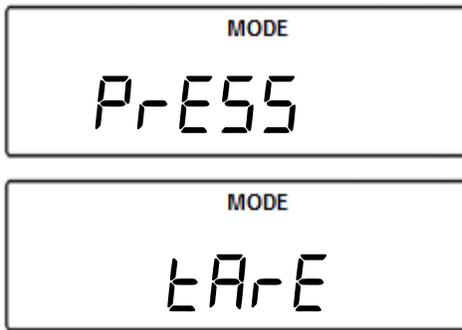

- Place an empty vial or counting container on the pan. The container’s weight will be displayed. Wait for the weight to stabilize and press the T key to tare.

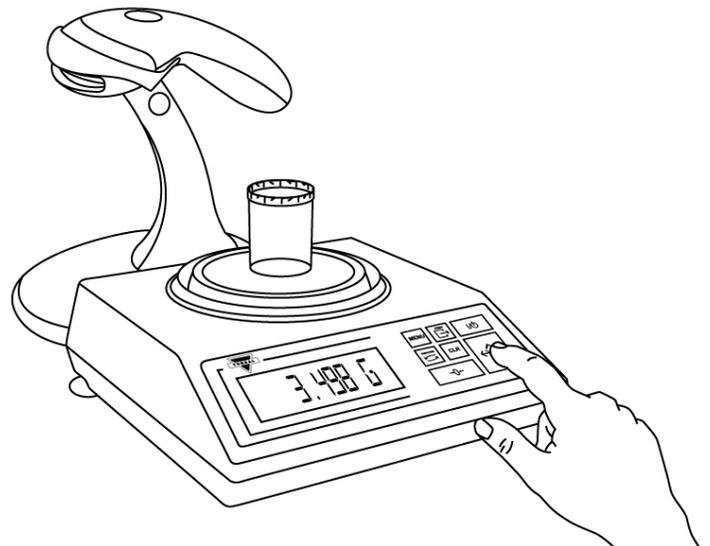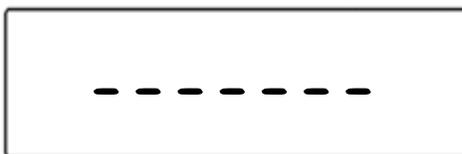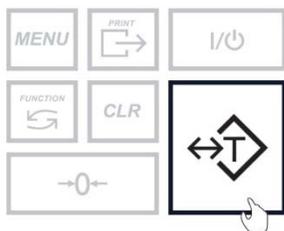


- Once the scale has finished taring “ScAn bottle” will appear on the display indicating you should scan the NDC code associated with the drug.

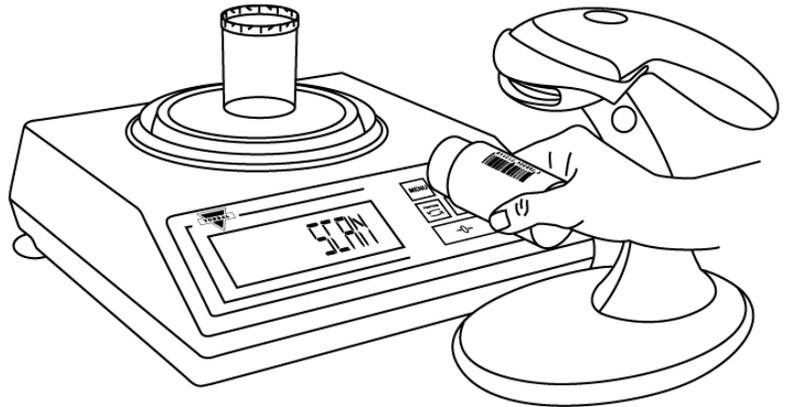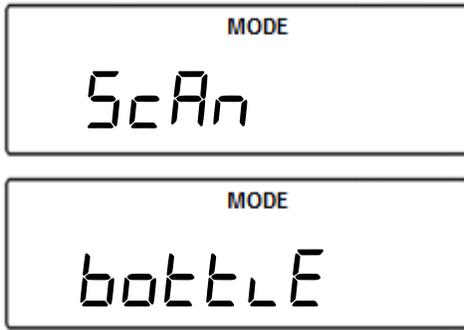

Note: If Rx Verification has been enabled, the bottle NDC and Rx NDC label must be scanned to allow the scale to check whether there is a match between the two codes.

- The scale will display FILL indicating to begin filling. Begin to pour the pills into the container.
- Once an accurate piece count has been taken, the container and it's content may be removed from the scale. After removing the container, done will appear on the screen.

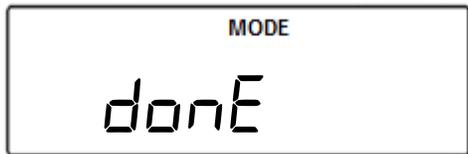

- To finish the count and begin a new one, press the Clear key.

