



Totalizing (TOTAL)

Totalizing is used to calculate the summation of multiple weights that are recorded sequentially. An “F” indicator will appear on the screen to indicate the scale is in Totalizing Mode.

1. To enter Totalizing Mode, press the “Function” key. Select “TOTAL” by pressing the “T” key.

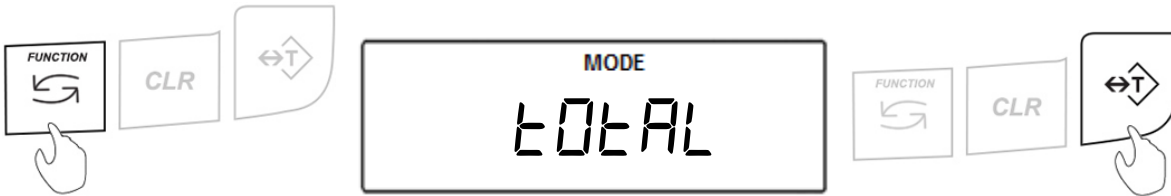


2. Place an empty weighing container on the pan. The container’s weight will be displayed. Wait for the weight to stabilize and press the “T” key to tare.

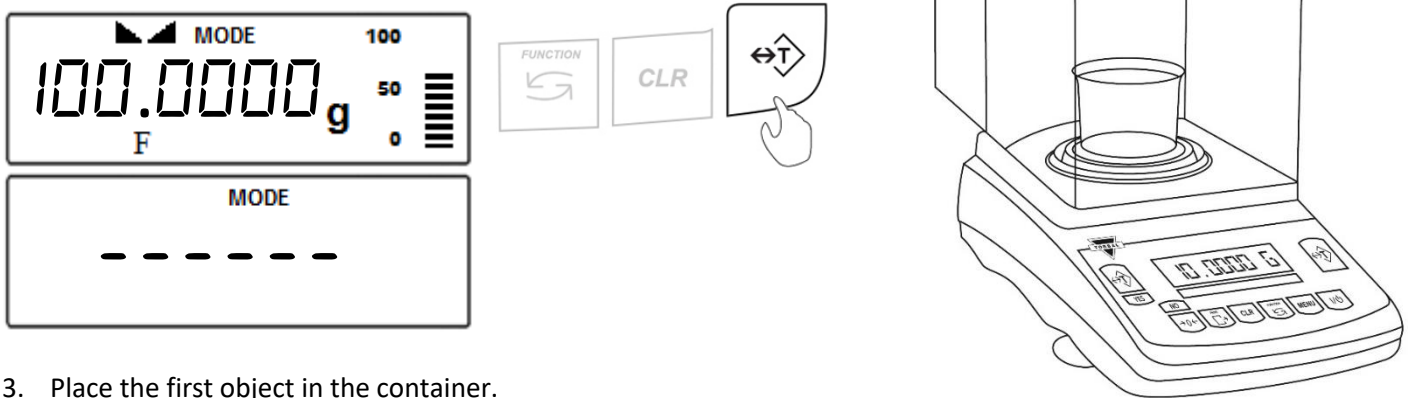

3. Place the first object in the container.


Automatic and Semi-Automatic:

When set to Auto or Semi-Auto, the weight of the object will be added to the total automatically. A TOTAL indicator will appear on the display.

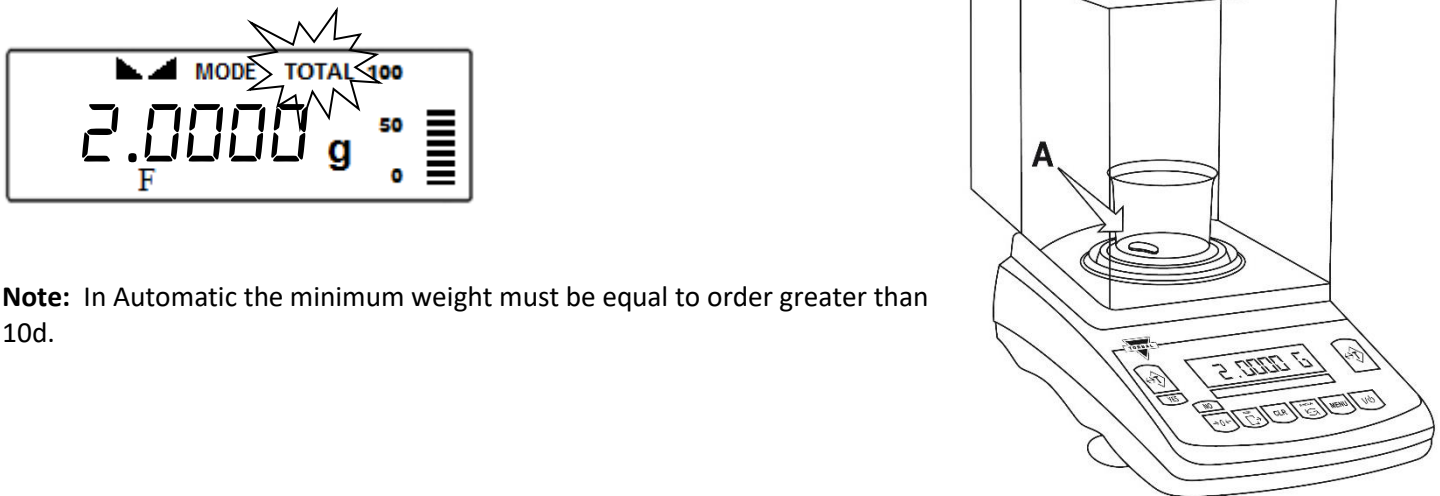

Note: In Automatic the minimum weight must be equal to or greater than 10d.

Manual: When set to Manual, press the “FUNCTION” key to add the weight to the total. A *TOTAL* indicator will appear on the display.

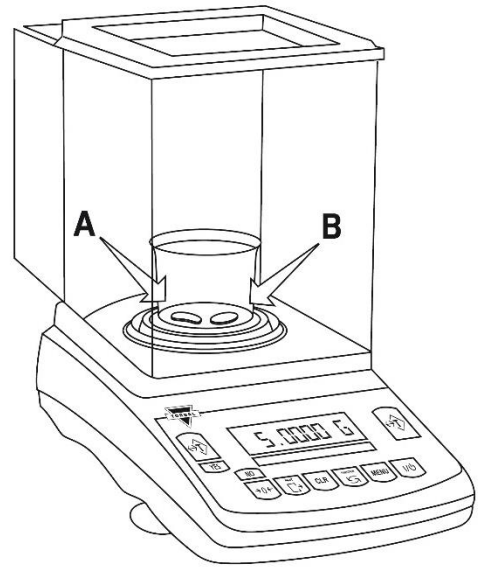


Note: With each added weight, the grand total result will appear on the display for approximately three seconds. To recall the grand total at any time, press the “MENU” key.

- Once a weight has been added to the total, remove the object from the pan or press the “T” key.



- Repeat **Steps 3 – 4** for the remainder of the samples.

- To clear the grand total and reset the function press the “CLR” key. “YES-NO” will appear on the display. To clear the function, press the “T” (Yes) key. To cancel the clear, press the “→0←” key.

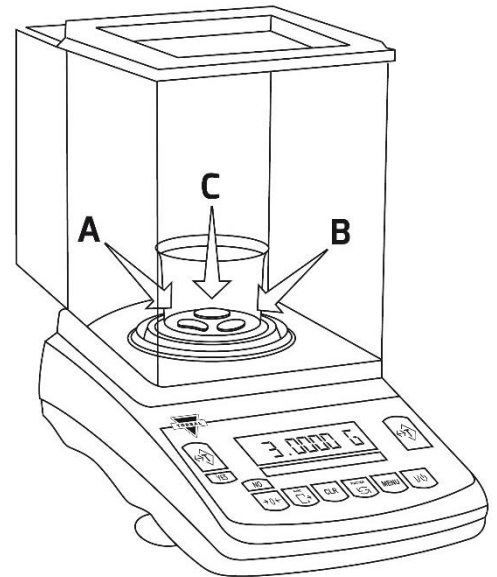

To ACCEPT

To CHANGE



- To obtain the total press the M key. To reset the compounding function and start a new recipe press the CLR key again while the total is displayed.

- To print the recipe, press the Print key.

