

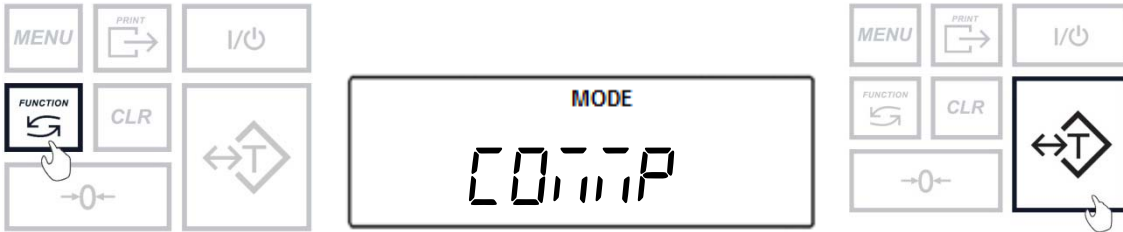


# Compounding


1. To enter Compounding Mode, press the Function key. Select “COMP” by pressing the T key.

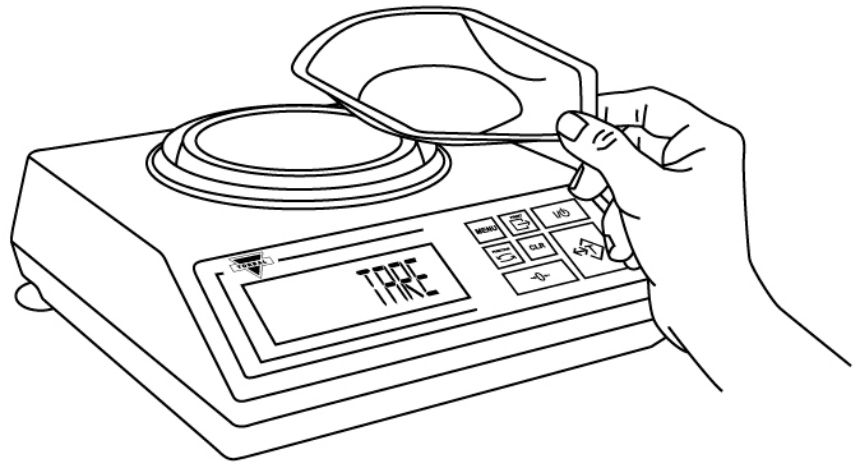



2. “PRESS TARE” will appear on the scale’s display, indicating to tare a container.

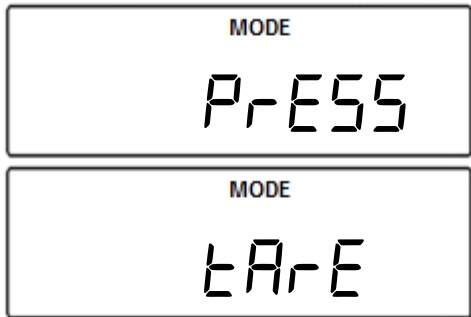


3. Place an empty container on the pan. The container’s weight will be displayed.

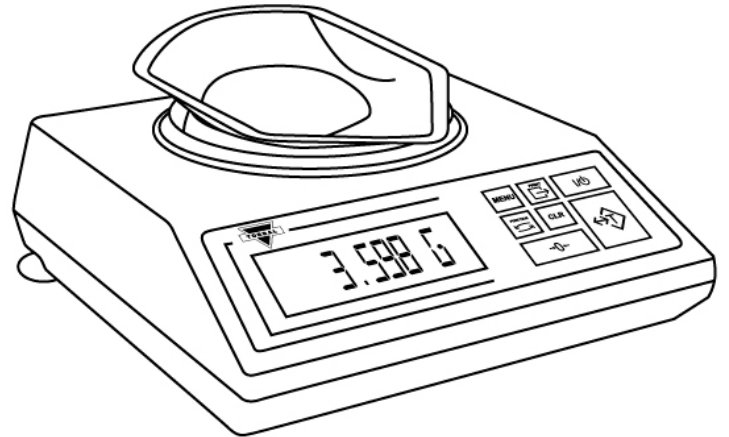


4. Wait for the weight to stabilize and press the T key to tare.

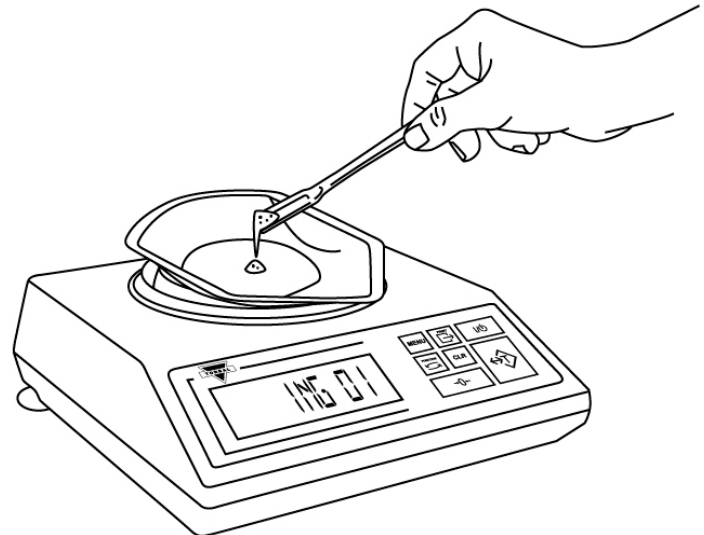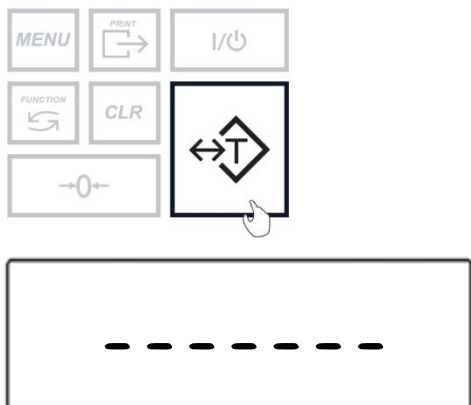



- After the container is tared, “1.9 01” will appear on the display, indicating the scale is ready for the first ingredient.

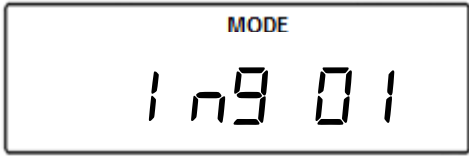



- Add the ingredient until the desired weight is displayed and the stability indicator appears on the display.

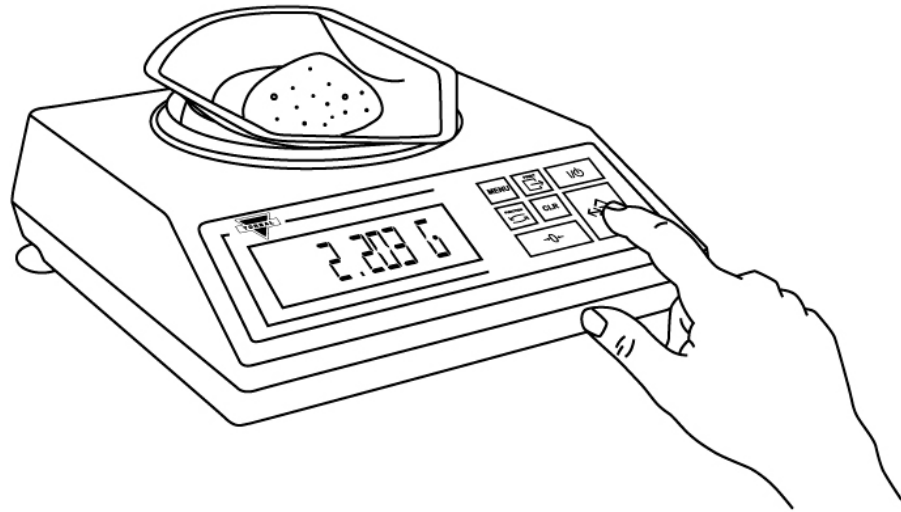

- Once a desired weight is displayed press the T key to store the ingredient in the scale’s memory.

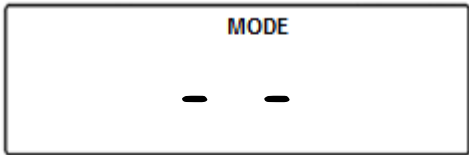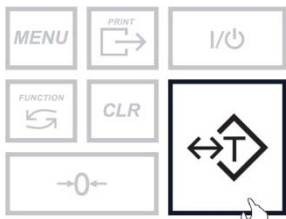

- Once the weight is stored, the scale will display “PRESS TARE” indicating to tare a container for the second ingredient. You may now repeat steps 3 to 7 to add more ingredients to the recipe.

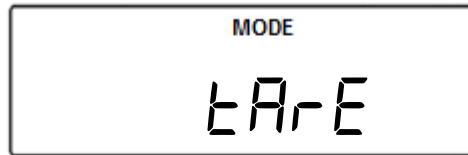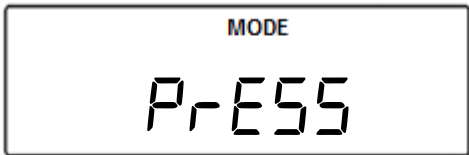

- To obtain the total press the CLR key. To continue adding more ingredients press the T key while the total is displayed. To reset the compounding function and start a new recipe press the CLR key again while the total is displayed.

