

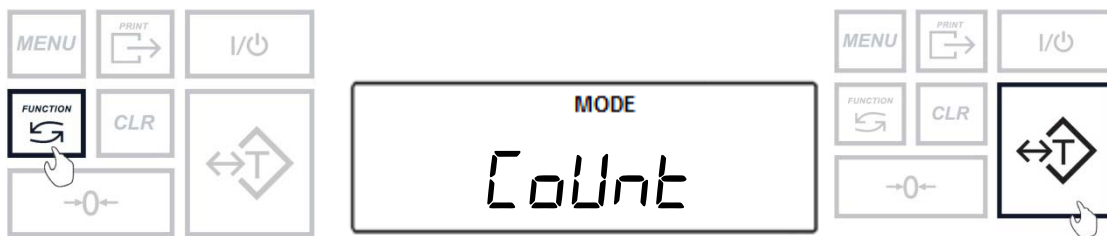


Pill Counting (DRX-4C2 Models)




1. To enter Pill Counting Mode, press the Function key. Select "Count" by pressing the T key.

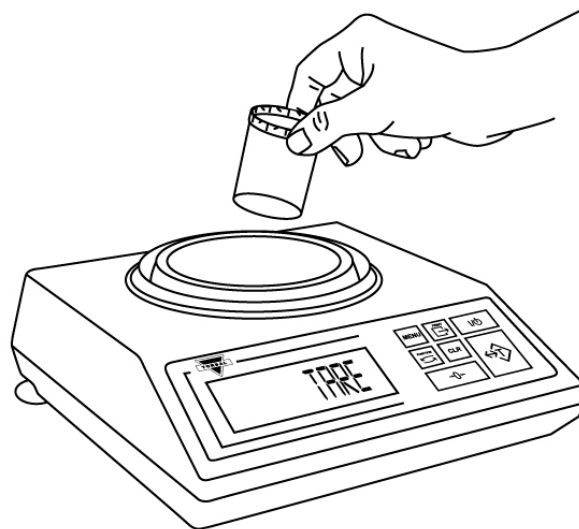



2. "PRESS TARE" will appear on the scale's display, indicating to tare a counting container.

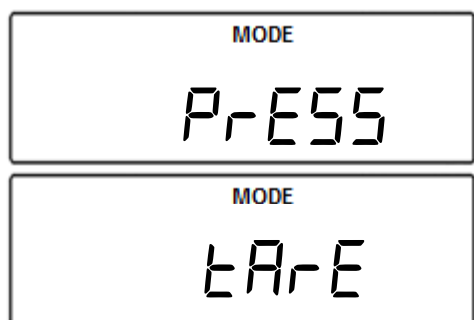




3. Place an empty vial or counting container on the pan. The container's weight will be displayed. Wait for the weight to stabilize and press the T key to tare.

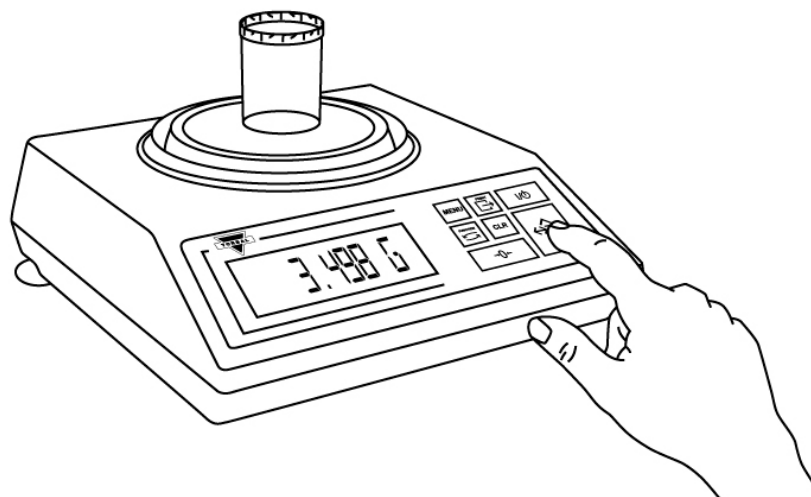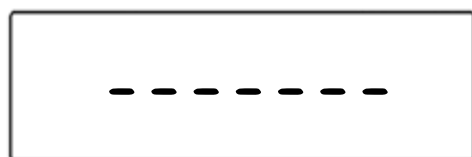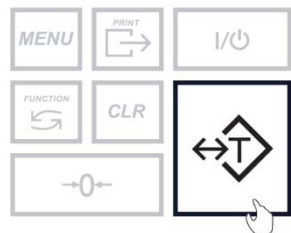



- The scale will display SAPL ON (SAPL ON). Manually count the required piece sample and place it in the container on the pan.

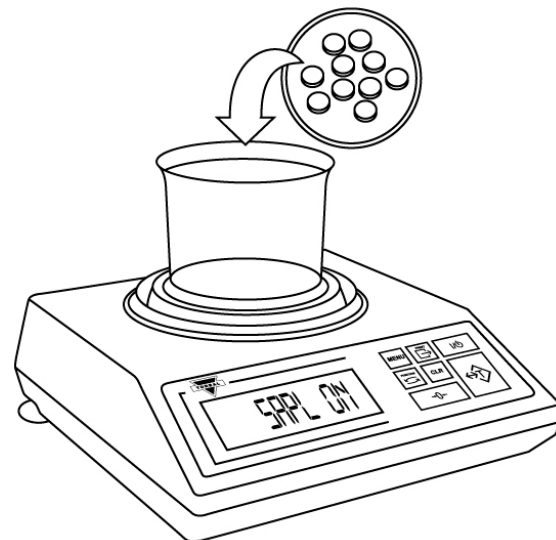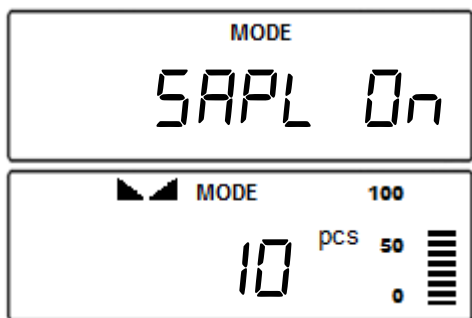

- The scale will display the weight of the sample. Press the T key to confirm the sample weight.

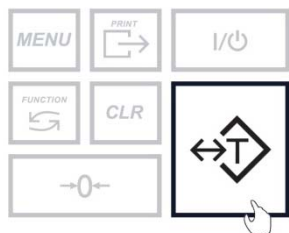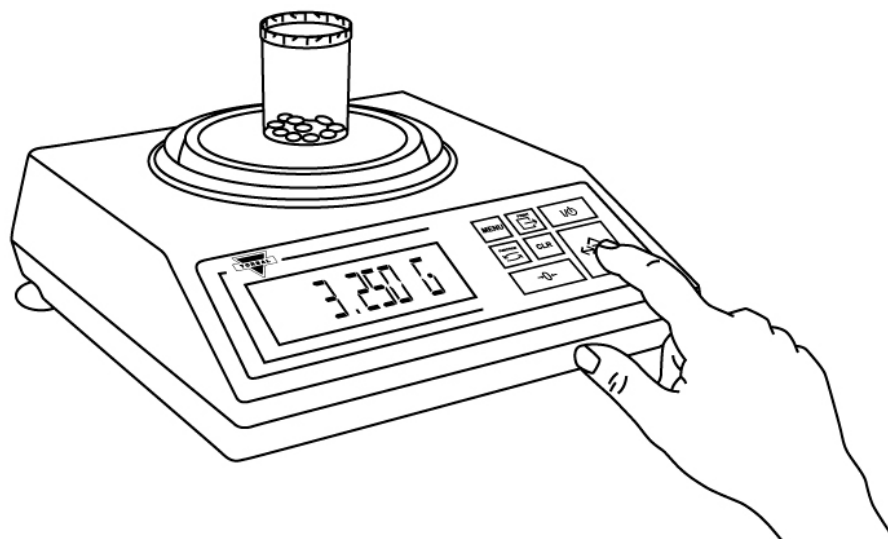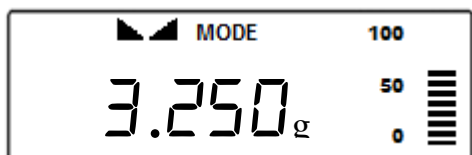

- The scale will display FILL indicating to begin filling the remainder of the count. Begin to pour the pills into the container.
- Once an accurate piece count has been taken, the container and its contents may be removed from the scale. After removing the container, *done* will appear on the screen.

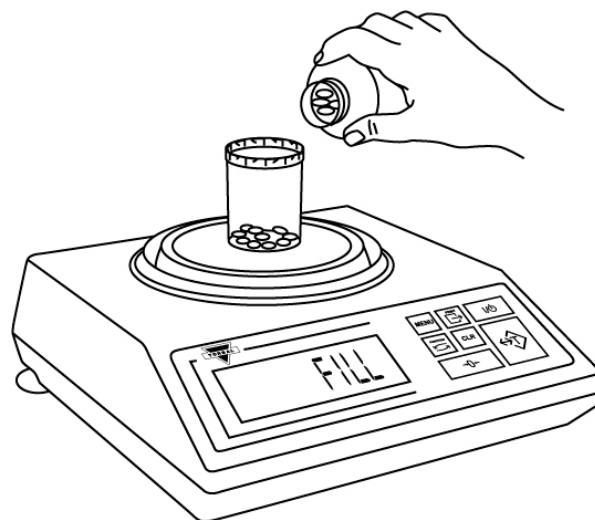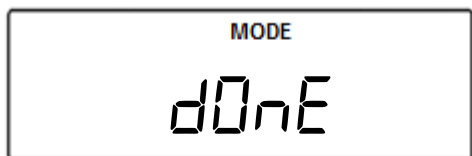

Note: If more pieces have been placed into the container than actually desired, you may use tweezers, a pharmacy spoon, or a similar instrument to remove the container from the pan, dispense out the excess pieces, and place it back on the pan to recheck the count

- To finish the count and begin a new one, press the Clear key.

