


# Animal Weighing

Before using this reference guide, enable and configure Animal Weighing by following the steps listed in Chapter 7.1.3 of the Instruction Manual. To enter the Animal Weighing mode press the “Function” key, and select the function by pressing the “T” key while “Animal” is displayed.

1. Place a container on the pan and tare it by pressing the “T” key.

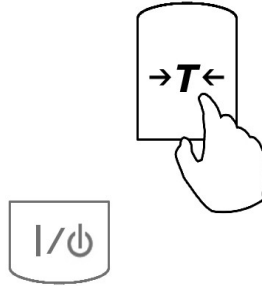

2. Place the animal into the container.




<i>Automatic and Semi Automatic modes</i>	<i>Manual mode</i>
Animal Weighing will begin <b>automatically</b> .	Press the "Function" key to initialize Animal Weighing.

3. The scale will count down a time interval.




4. After the time interval has elapsed the result will be displayed.

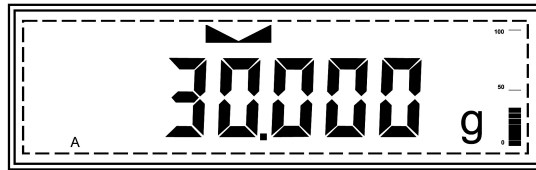


5. The animal can be removed from the pan



*Automatic mode:* Result will be automatically cleared; the scale will perform an automatic tare.

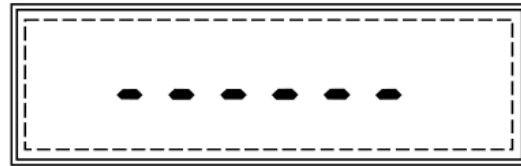

*Manual and Semi Automatic modes:* To clear the result, press the “C” key. After clearing the function press “T” to tare.

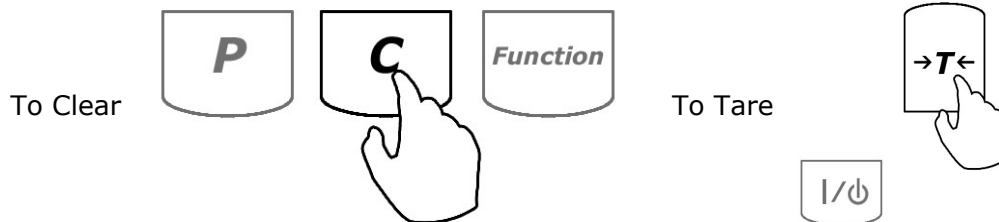

**For a detailed description of this feature please turn to Chapter 8(A) of the instruction manual.**

For technical questions or assistance please call toll free: 1-866-473-6900

Fulcrum Inc.  
 100 Delawanna Ave., Suite 120  
 Clifton, NJ 07014  
 (973) 473-6900  
<http://www.torbal.com>