



Display Hold


Before using this reference guide, enable and configure Display Hold by following the steps listed in Chapter 7.1.3 of the Instruction Manual. To enter the Display Hold mode press the “Function” key, and select the function by pressing the “T” key while “Display” is displayed.

1. Place a container on the pan and tare it by pressing the “T” key.

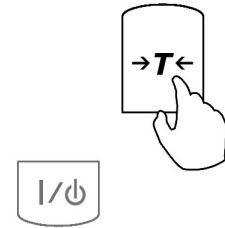

2. Place the weight in the container located on the pan.

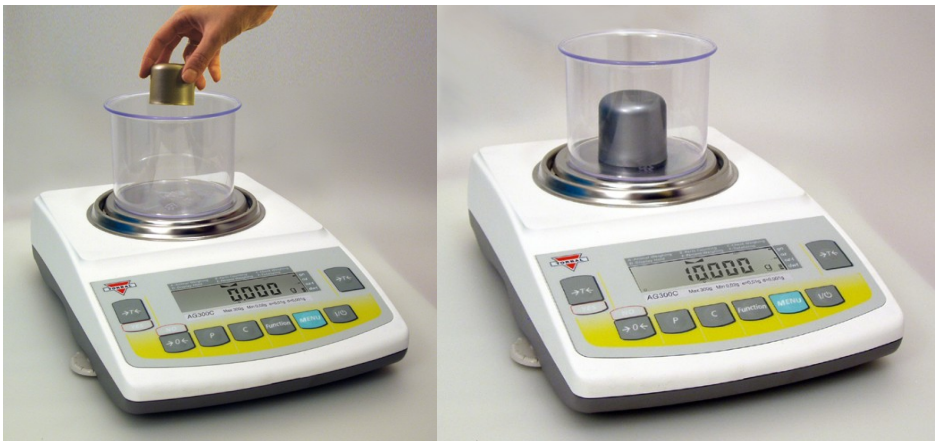



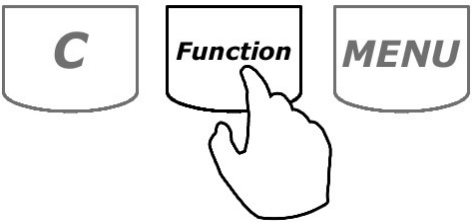
Automatic and Semi-Automatic	Manual
<p>The scale will automatically lock the weighing result on the display as soon as it stabilizes.</p>	<p>Display hold is initialized by pressing the “Function” key. Once the weight is stable press “Function” to lock the result on the display.</p> 



AGC Series: Quick Reference Guide

3. Remove the weight from the pan.

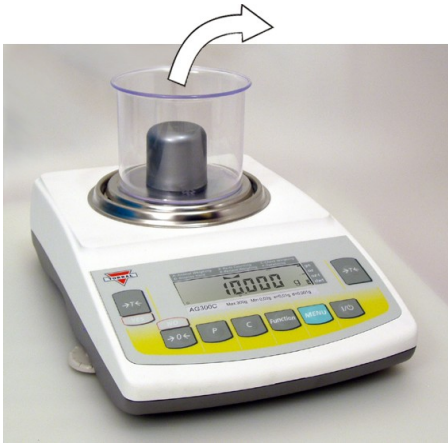

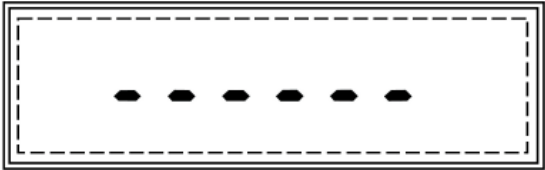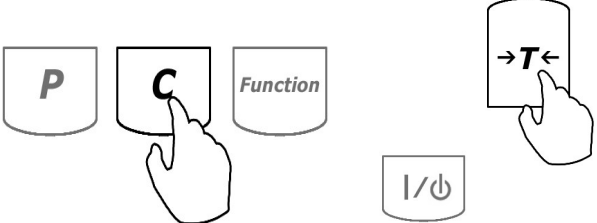Automatic	Semi-Automatic and Manual
<p>The result will be automatically cleared after the weight is removed from the pan. The scale will perform an automatic tare and prepare for the next weighing.</p> 	<p>To clear the weighing result press the “C” key. After the result has been cleared the scale must be tared. To tare the scale press the “T” key.</p> 

For a detailed description of this feature please turn to Chapter 8(D) of the instruction manual.

For technical questions or assistance please call toll free: 1-866-473-6900

Fulcrum Inc.
100 Delawanna Ave., Suite 120
Clifton, NJ 07014
(973) 473-6900
<http://www.torbal.com>