




Parts Counting

Before using this reference guide, enable and configure Parts Counting by following the steps listed in Chapter 7.1.3 of the Instruction Manual. To enter the Parts Counting mode press the “Function” key, and select the function by pressing the “T” key while “Parts” is displayed.

1. Place a container on the pan and tare it by pressing the “T” key.

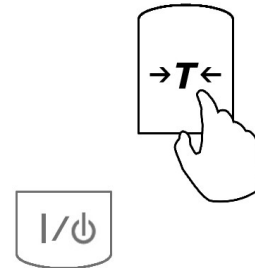



2. Confirm the sample size by pressing the “T (Yes)” key.

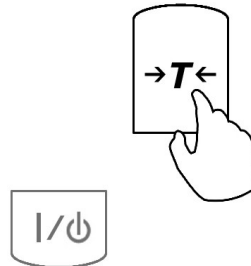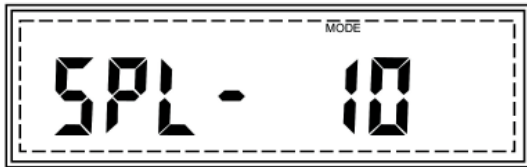


3. Manually count out the sample and place it in the container located on the pan.



4. Press the “T (YES)” key to confirm and accept the sample weight.

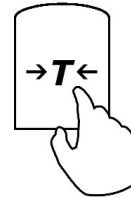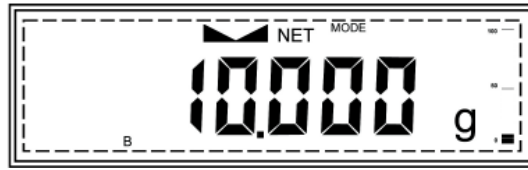

5. Begin to fill the remainder of the pieces.

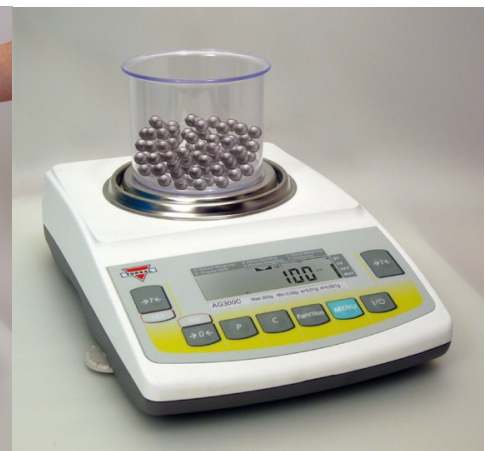

For a detailed description of this feature please turn to Chapter 8(B) of the instruction manual.

For technical questions or assistance please call toll free: 1-866-473-6900

Fulcrum Inc.
100 Delawanna Ave., Suite 120
Clifton, NJ 07014
(973) 473-6900
<http://www.torb.com>