


# Totalizing



Before using this reference guide, enable and configure Totalizing by following the steps listed in Chapter 7.1.3 of the Instruction Manual. To enter the Totalizing mode press the “Function” key, and select the function by pressing the “T” key while “Totalize” is displayed.

1. Place a container on the pan and tare it by pressing the “T” key.

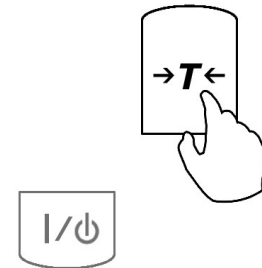


2. Place the first object in the container located on the pan.

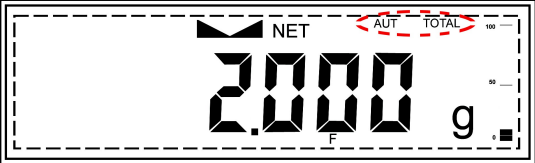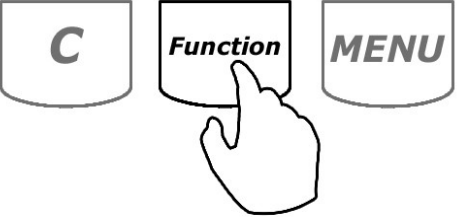


Automatic	Manual
<p>The weight of the object will be added automatically. Once the weight stabilizes, the <b>TOTAL</b> indicator will appear on the display, signaling that the weight has been added to the grand total</p> 	<p>Once the weight stabilizes and the stability indicator appears on the display press the “Function” key to add the weight to the grand total.</p>  <p>Add to Total</p>


With each added weight, the grand total result appears on the display for approximately three seconds. To recall the grand total at any time, press the “Menu” key.

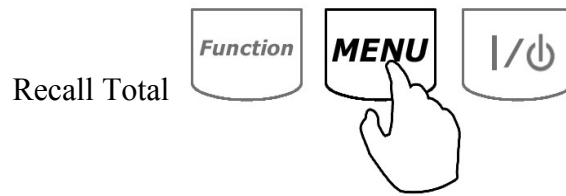


- Once a weight of an object has been added to the total, remove the object from the pan or press “T” to tare the scale.

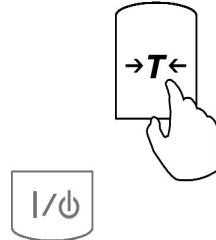

- After the scale has been tared, Steps 1 through 3 can be repeated in order to add more weighing results to the grand total.

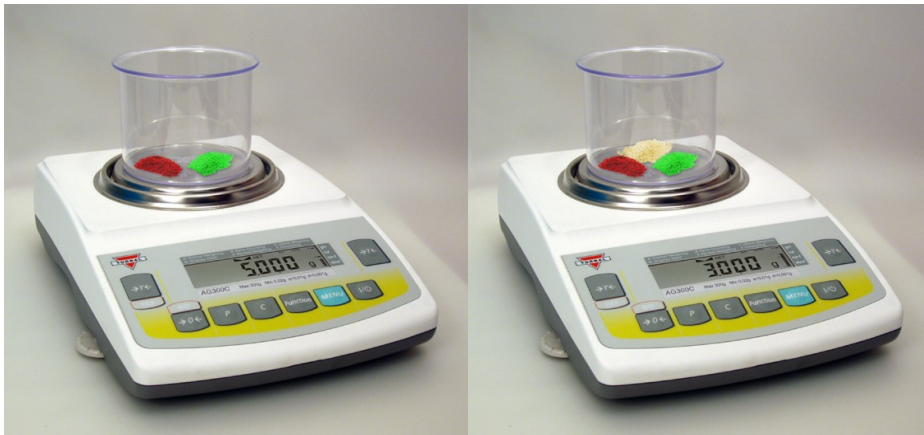

- To clear the grand total and reset the function press the “C” key.

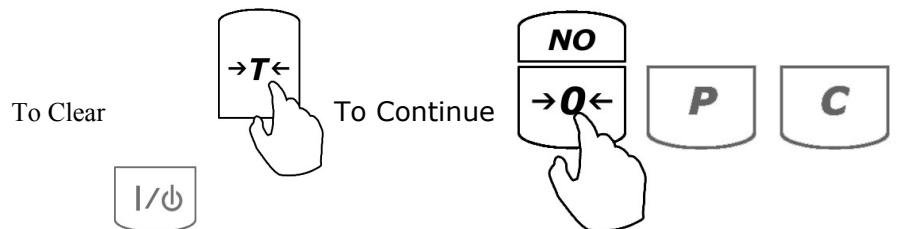

**For a detailed description of this feature please turn to Chapter 8(F) of the instruction manual.**

For technical questions or assistance please call toll free: 1-866-473-6900

Fulcrum Inc.  
 100 Delawanna Ave., Suite 120  
 Clifton, NJ 07014  
 (973) 473-6900  
<http://www.torbald.com>