

Pan Installation Guide

Models: ATS & BTS



1. Carefully remove the Moisture Analyzer and all its components from the packaging and place them on a stable base where the unit will not be affected by any mechanical vibrations or air movements.
2. After removing the Pan Shield (4), Pan Support (5), and Pan Handle (6) from their packaging, open the Drying Chamber (3) and carefully install the Pan Shield (4) on the three Pan Shield Posts (8).
3. Once the Pan Shield is installed on the shield posts, gently insert the Pan Support (5) into the opening of the Analyzer's mechanism.
4. After installing the Pan Support, place a Disposable Pan (7) on the Pan Handle (6), and then place it on the Pan Support so that the handle rests on the Pan Shield without touching the Pan Support.

