

## Packing Specifications


Please retain the original packaging and use it whenever storing or transporting the analyzer. Always pack the analyzer as shown below. When packing the unit, be sure to remove the pan, pan base, draft ring and the chamber base. Shipping the analyzer with the pan components installed may damage the unit. Always place the items in their designated compartments, as shown below.

### Pan Components

Pan Shield
Pan Support
Pan Handle
Disposable Pans
Sodium Tartrate Dihydrate
Forceps



### Scale

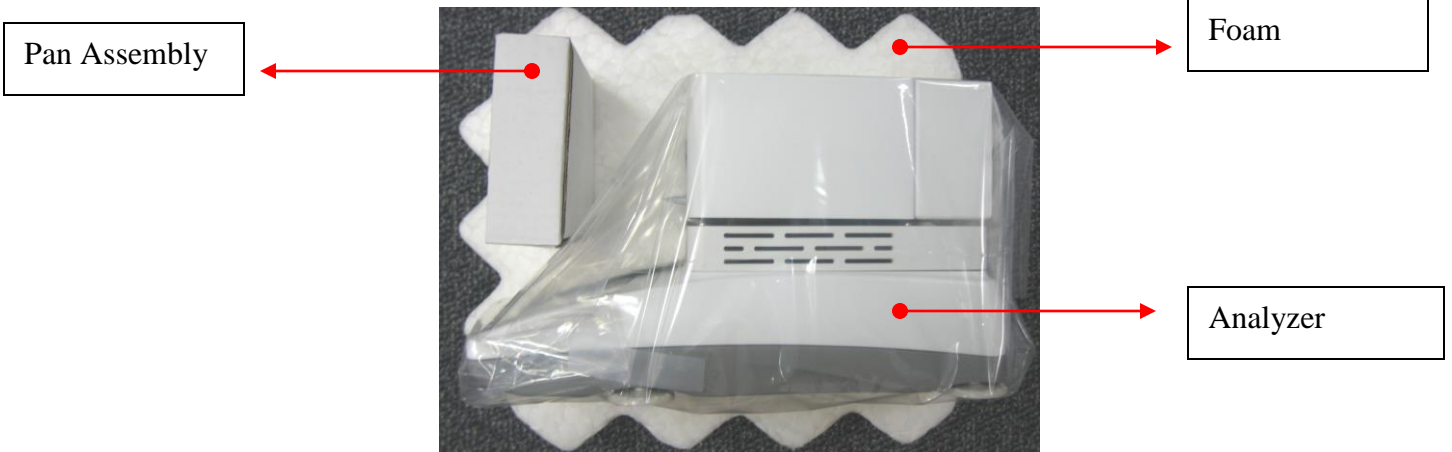

**- Never Ship or Transport the Analyzer with Pan Components Inside of the Drying Chamber! -**