
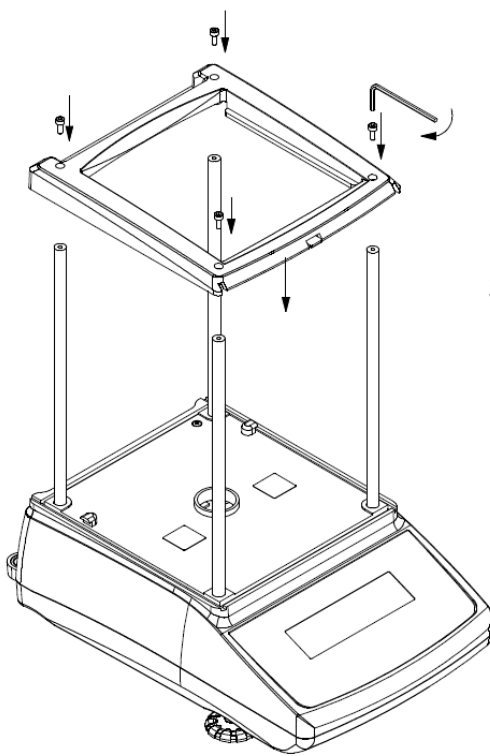


Getting Started


ATN60A | AT110A


1. Carefully remove the scale, pan, and all components out of the packaging. Place them on a stable surface where the scale will not be compromised by any mechanical vibrations or high air movements.
2. Firmly screw in the chamber posts into the screw hole openings located in the corners of the scale's top housing.

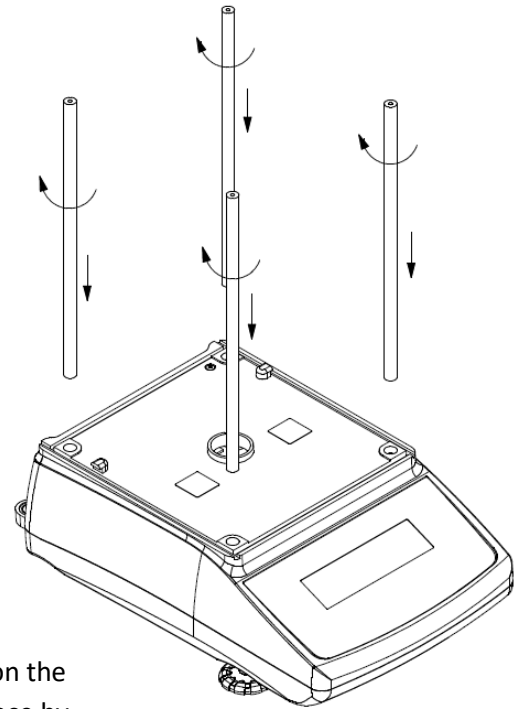

3. Place the chamber top frame on the posts and firmly secure it in place by screwing in the post crews with the provided wrench.



4. Install the back-chamber glass by inserting the bottom edge into the scale housing (C') and securing the top edge in the chamber top frame (C'').
5. Install the left and right chamber access door by sliding it into the scale housing and the chamber top frame (C''').

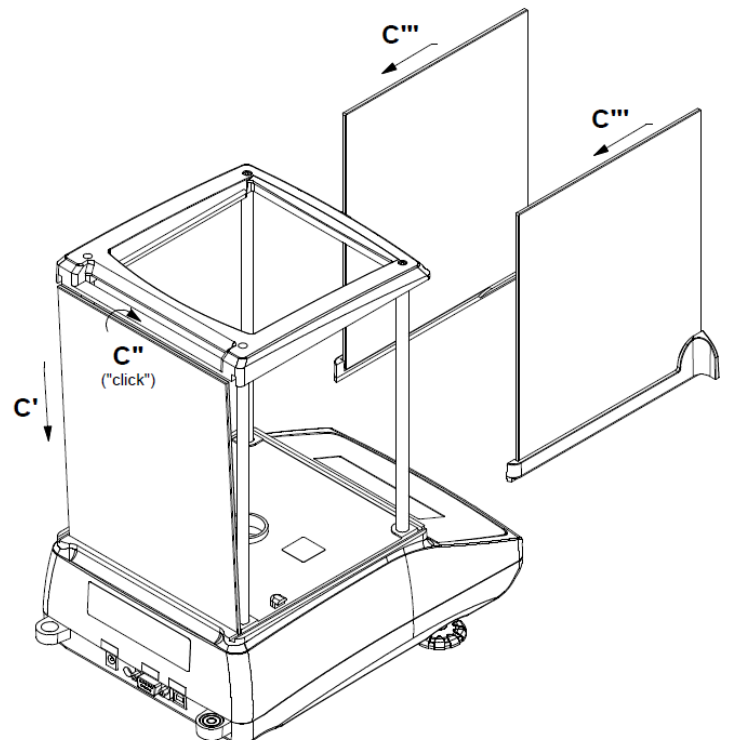

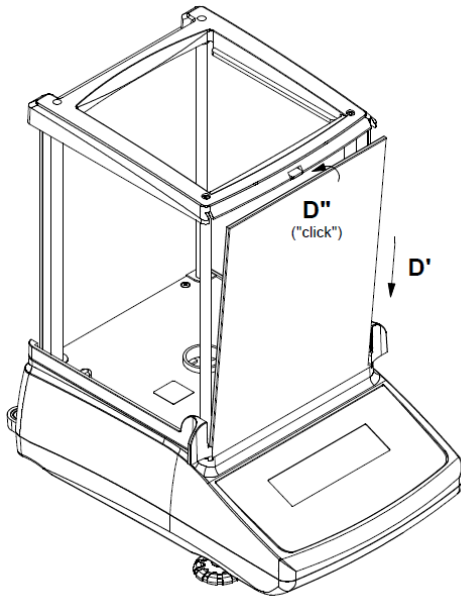

6. Install the front-chamber glass by inserting the bottom edge into the scale housing (D') and securing the top edge in the chamber top frame (D'').

7. Insert the top chamber glass access door by sliding in the rear edge of the glass onto the chamber top frame. (E')

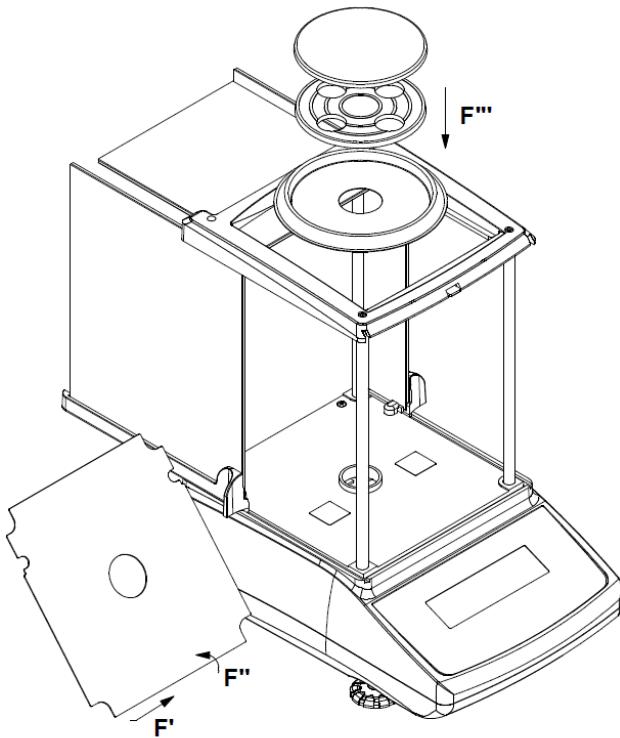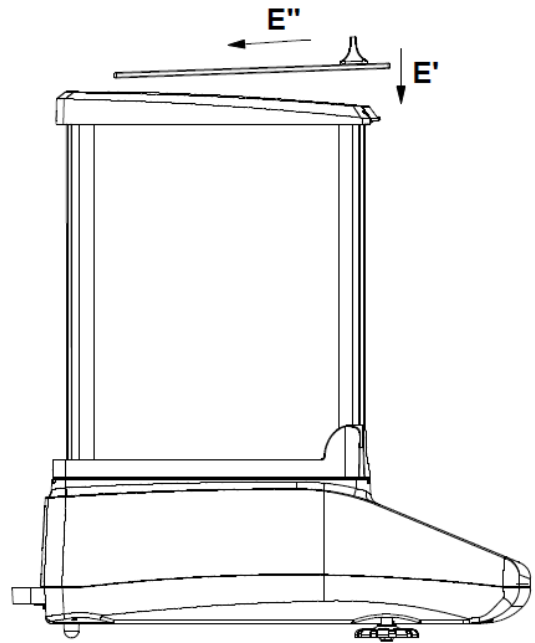

8. Install the chamber base (F') on the top of the scale's housing.
9. Once the chamber bases in place, install the draft ring, pan support and the pan (F''').