




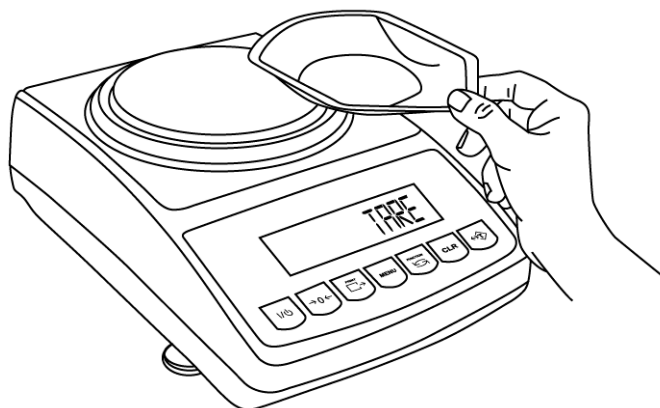


Compounding



1. "PrESS tArE" will appear on the scale's display, indicating to tare a container.

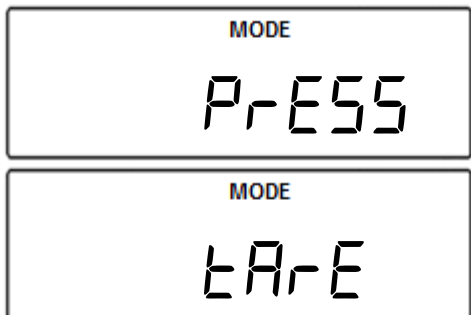


2. Place an empty container on the pan. The container's weight will be displayed.

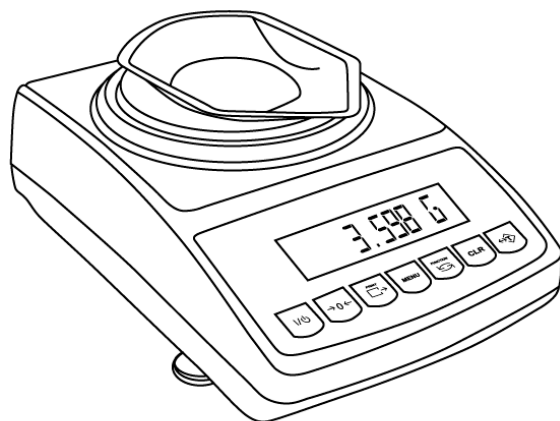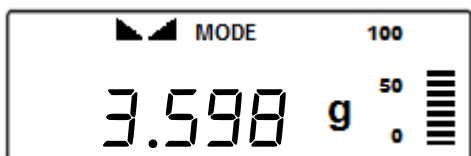


3. Wait for the weight to stabilize and press the T key to tare.

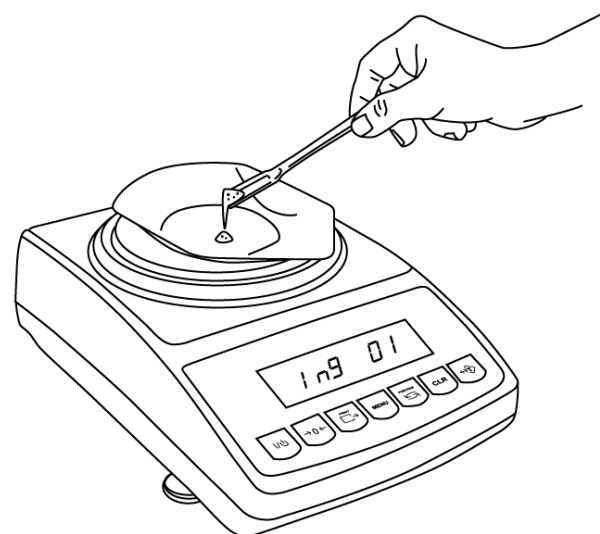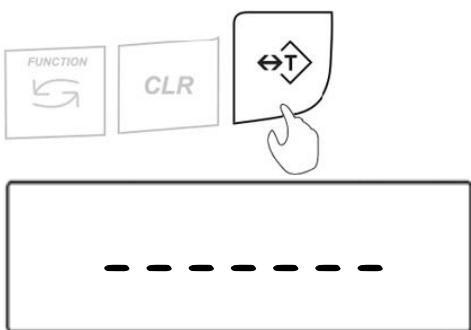


4. After the container is tared, "1.9 01" will appear on the display, indicating the scale is ready for the first ingredient.

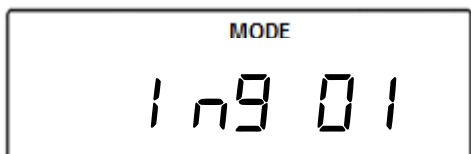



5. Add the ingredient until the desired weight is displayed and the stability indicator appears on the display.

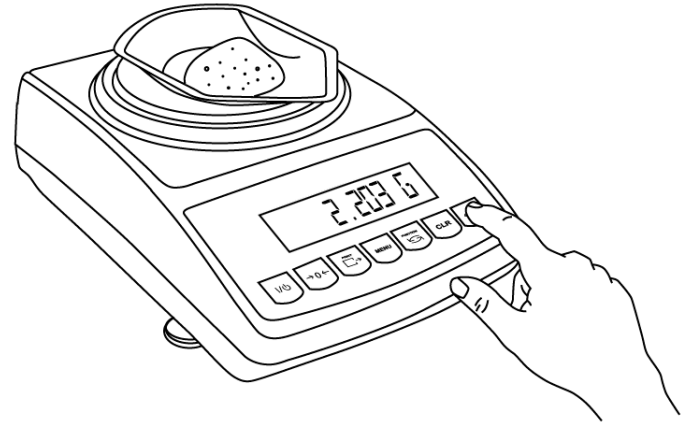


6. Once a desired weight is displayed press the T key to store the ingredient in the scale's memory.

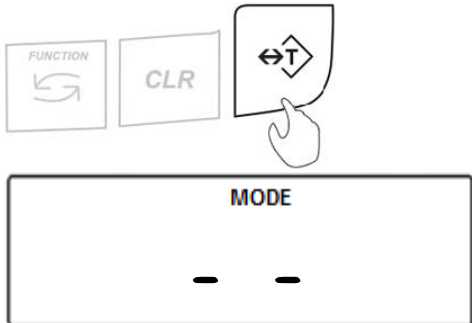

7. Once the weight is stored, the scale will display "P-R-E-S-S T-A-R-E" indicating to tare a container for the second ingredient. You may now repeat steps 3 to 7 to add more ingredients to the recipe.

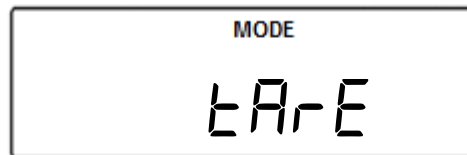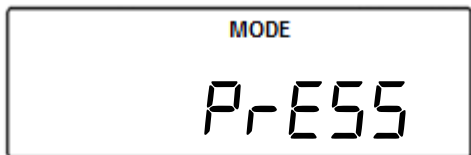

8. To obtain the total press the CLR key. To continue adding more ingredients press the T key while the total is displayed. To reset the compounding function and start a new recipe press the CLR key again while the total is displayed.

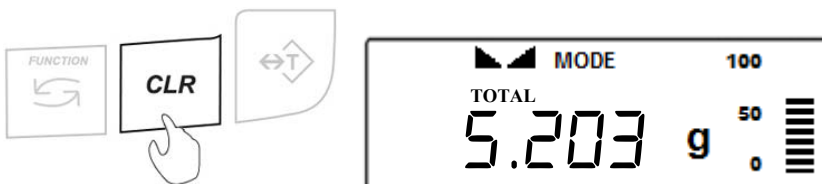

9. To print the recipe, press the Print key.

